



### □ G1540CVR Portable Eyewash Cover

## Assembly Instructions

1. Fully empty unit. Allow all water to dispense through dispenser before attempting to lift. Use 2 persons to lift unit when emptying.
2. Open the cover and spread out as shown in Figure 1.
3. Place eyewash unit into cover as shown in Figure 2. Pull front of cover up and over the front "lip" of the tank.
4. Pull sides up to cover unit, keeping the center panel separated. With both sides elevated, secure Velcro strap on top of unit.
5. Situate the eyewash unit to best fit inside the cover and adjust the Velcro strap to obtain a secure fit. See Figure 3.
6. With activating panel of eyewash unit in the down position, attach approximately 1 inch of the Velcro from the pull down flap to the Velcro on the activating panel as picture in Figure 4 and Figure 5.
7. Lift activation panel of eyewash unit to the "ready" or "up" position with pull down flap attached. See Figure 6.
8. Starting from the bottom and working up, secure Velcro along the edges of the pull down flap on the cover. The unit is now protected.
9. Fill unit as per the original instruction provided with the eyewash unit.

## Operating Instructions

The unit is operated in one easy step. Simply pull down on the dust cover handle (located at the top of the yellow activating panel) in one swift motion. The dust cover flap will fall away and the yellow activating panel will rest in the basin of the unit initiating the eyewash flow.

## Cleaning Instructions

The dust cover may be cleaned with mild soap and water.

Distributed by:  
Safety Emporium  
PO Box 1003  
Blackwood, NJ 08012

Ph: (866) 326-5412 toll-free  
Fax: (856) 553-6154  
esupport@safetyemporium.com  
www.safetyemporium.com

## Installation Photos

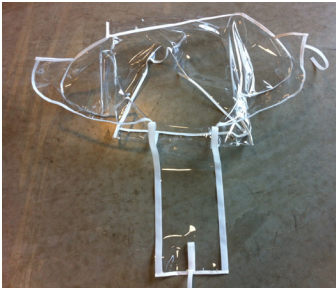


Figure 1



Figure 2



Figure 3



Figure 4



Figure 5



Figure 6



Complete - Front View



Complete - Top View